



**QAQC LAB**

www.qclabequipment.com



## Powder Trier

High Quality Trier - ideal for powders, granules & grains

The QAQC LAB Powder Trier is ideal for taking a large volume, cross sectional sample from a container. The sampler has a robust all stainless steel construction and is suitable for sampling cereals, grains, free flowing powders, granules and even slightly cohesive powders

- Cross section (composite sample)
- 316 stainless steel, bright finish
- Suitable for taking a large volume sample from compacted powders, grains etc.

The cutting scroll helps sampler go into product

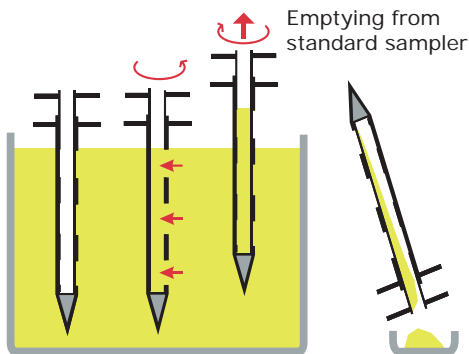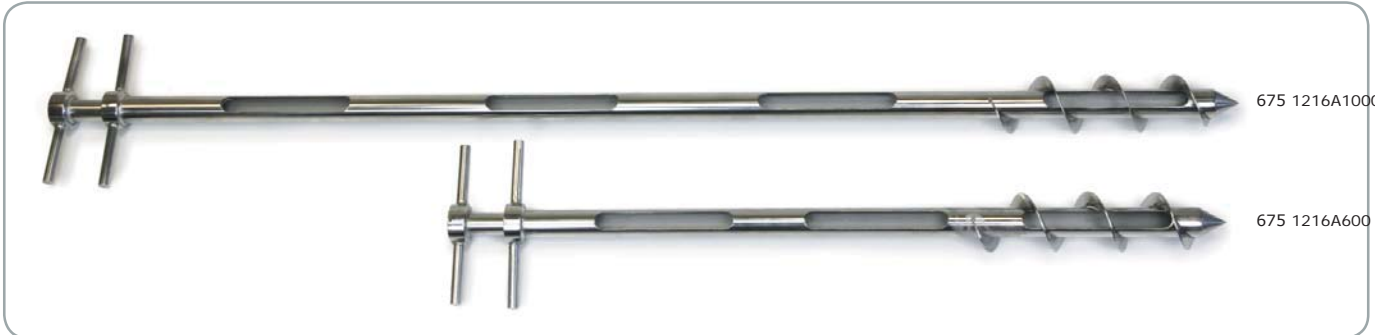

### Operation

1. Insert the sampler into the product
2. Turn the handle to open the slot - the product can now be collected in the central tube
3. Turn the handle to close the slot
4. Remove the sampler
5. Invert the sampler to discharge the sample

Note: The cutting scroll is not shown in the sketches

### QUICK QUOTE

### ONLINE PRICING

Stock No.	Length	Number of Slots	Max Sample Volume
675 1216A600	600mm	3	110ml
675 1216A1000	1000mm	4	190ml
675 1216A1500	1500mm	5	300ml