



QAQC LAB


www.qclabequipment.com

U-D Cohesive Sampler

Unit Dose Sampler for Cohesive Powders


A popular sampler in the pharmaceutical industry for taking low volume (unit dose) cohesive powder samples. Allows accurate samples to be taken.

- Single Point Unit Dose
- 316 stainless steel
- Suitable for use with sticky, cohesive powders

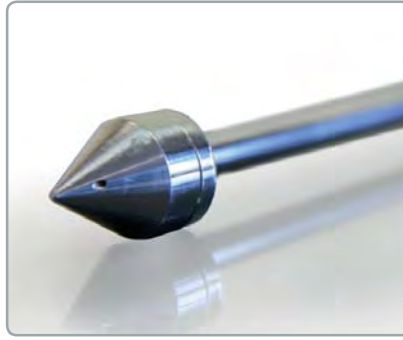

Operation

1. Insert the sampler into the product.
2. At the required depth push down the handle and rotate it through 180°.
3. Pull up sampler slightly (to fill the cell), then rotate handle 180° (to lock cell in closed position).
4. Remove sampler

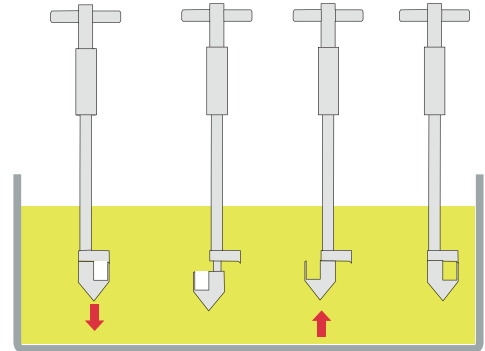

Sample Collecting Cells are available in a range of standard sizes.



25ml



50ml



75ml



100ml

QUICK QUOTE

ONLINE PRICING

Stock No.	Description
675 2000S01	U-D Cohesive Sampler Set - consisting of: 1 off Handle, Sample Cells - 1 each of: 0.25ml, 0.5ml, 0.75ml and 1.0ml 1 off Standard Sampler Case
675 2000A1000	Sampler body and handle - 1000mm length
675 2000T025	Sample Collecting Cell - 0.25ml
675 2000T050	Sample Collecting Cell - 0.5ml
675 2000T075	Sample Collecting Cell - 0.75ml
675 2000T100	Sample Collecting Cell - 1.0ml
675 8020N1100	Cleaning Brush - 1100mm length - narrow style

